

REES SCIENTIFIC Quick Reference

Responding to Centron Alarms

When called...

1. Listen for Alarm details.
2. Press 0 to acknowledge alarm.
3. Enter telephone ID code and #.
4. Listen & follow voice prompts to inhibit Alarm.
5. Press 1 for Global status
 - 2 to Check single input status
 - 3 to Inhibit an input
 - 4 to Enable an input
 - 5 to End Call
6. Listen & follow voice prompts for appropriate action.
7. Press 5 to End Call.

See Reverse side for Calling in to System.

Always

Press # after digits for system to accept entry.
End call by pressing 5. Do NOT just hang up.
Report activity to your supervisor & document actions.

For complete instructions visit www.reesscientific.com
or call Tech Support at 609.406.0073

REES SCIENTIFIC Quick Reference

Responding to Centron Alarms

When calling in...

1. Call System

Node	Phone #
Node	Phone #
Node	Phone #
Node	Phone #
Node	Phone #
Node	Phone #
Node	Phone #
Node	Phone #
Node	Phone #
Node	Phone #
Node	Phone #
Node	Phone #
2. Enter telephone ID code and #.
3. Press 1 for Global status
 - 2 to Check single input status
 - 3 to Inhibit an input
 - 4 to Enable an input
 - 5 to End Call
4. Listen & follow voice prompts for appropriate action.
5. Press 5 to End Call.

See Reverse side for when System calls.

For complete instructions visit www.reesscientific.com
or call Tech Support at 609.406.0073